

Biorbyt dNTP-Mutagenesis Kit

Cat#: orb533432 (User manual)

Biorbyt dNTP-Mutagenesis Kit Random Mutagenesis by dNTP Analogs

For in vitro use only! Shipping: shipped on gel packs Storage Conditions: store at -20 °C Additional Storage Conditions: avoid freeze/thaw cycles, aliquoting of the modified nucleotides prior to use is strongly recommended Shelf Life: 12 months

Description:

Biorbyt Mutagenesis Series

Within three billion years of evolution, nature has produced a plethora of proteins simply by repeated cycles of random mutagenesis followed by *in vivo* selection for superior function of the encoded proteins. This example of natural evolution has guided researchers within the last two decades to develop strategies for *in vitro* permutation of proteins.

Among the variety of strategies applied, three major powerful techniques have emerged.

Random Mutagenesis by dNTP Analogs

This method is based on the incorporation of mutagenic dNTP analogs, such as 8-oxo-dGTP and dPTP, into an amplified DNA fragment by PCR. The mutagenic dNTPs are eliminated by a second PCR step in the presence of the four natural dNTPs only, resulting in a rate of mutagenesis of up to 20%.

→Biorbyt dNTP-Mutagenesis

Random Mutagenesis by Error-Prone PCR

Developed by Caldwell & Joyce (1992) this method introduces mutations in the gene of interest using a PCR reaction under conditions that induce an increased error-rate of the DNA-polymerase. The rate of mutagenesis achieved by error-prone PCR is in the range of 0.6-2.0%.

→Biorbyt Error-Prone Kit

Random Mutagenesis by DNA Shuffling

Developed by Stemmer (1994) DNA shuffling generates libraries by random fragmentation of one gene or a pool of related genes, followed by the reassembly of the fragments in a self-priming PCR reaction. This method allows the recombination of sequences from different, related genes. The overall rate of mutagenesis is approx. 0.7%.



\rightarrow Biorbyt DNA-Shuffling Kit

Jena Bioscience now offers all components necessary for each of these techniques 'ready-to-go' in a separate kit, accompanied by a streamlined documentation that maximizes success.

Content:

Taq Polymerase (red cap) 5 units/μl, 40 μl Mutagenesis Buffer (green cap) 10x concentration, 200 μl dNTP Mix (white cap) 10 mM each dNTP (dATP, dCTP, dGTP,dTTP), 100 μl dPTP (yellow cap) 10 mM, 40 μl 8-oxo-dGTP (blue cap) 10 mM, 40 μl PCR-grade Water (white cap) 2x 1 ml

Random mutagenesis by dNTP analogs

8-oxo-dGTP as well as dPTP are mutagenic dNTP analogs which are incorporated into DNA by PCR using Taq polymerase [Zaccolo *et al.*, Zang *et al.*].

8-Oxo-dGTP is incorporated opposite template adenine yielding two transition mutations: $A \rightarrow C:T \rightarrow G$. A total rate of mutagenesis of approx. 2% (Fig. 2) has been reported, with $A \rightarrow C:T \rightarrow G$ occurring at a ratio of approx. 1:1.5.

The dNTP analog dPTP has been reported to be approx. 10-fold more mutagenic than 8-oxo-dGTP (Fig. 2). The mutations induced by dPTP occur at a ratio of approximately 5:4:1:1 ($A \rightarrow G:T \rightarrow C:G \rightarrow A:C \rightarrow T$), with a total rate of mutagenesis of up to 19%. Random mutagenesis induced by 8-Oxo-dGTP and/or dPTP is carried out in a two-step PCR process. First, the target DNA fragment is amplified in the presence of the four natural dNTPs plus the mutagenic analogs dPTP and/or 8-Oxo-dGTP. The rate of mutagenesis can be easily controlled by the number of PCR-cycles (Fig. 2).

The product of the first PCR is then subjected to a second PCR in the absence of mutagenic analogs. This step eliminates the non-natural analogs from the target DNA before cloning and transformation.

Recommended assay preparation

- For a 50 μ l reaction, take 5 μ l of 10x Mutagenesis Buffer in a sterile vial
- Add up to 25 fmol template DNA and 0.2-1 μM of appropriate primers
- Add 2.5 μl of dNTP Mix
- Add 2.5 µl of each mutagenic analog
- Add 1 µl (5 u) of Taq Polymerase
- Add PCR-grade Water to a final volume of 50 μl



Recommended thermo cycling conditions

Denaturation 92°C 1 min Annealing 55°C 1.5 min Extension 72°C 5 min Number of cycles: 5-30 (refer to Fig. 2) The rate of mutagenesis mainly depends on the number of cycles (Fig. 2) and can additionally be fine-tuned by the amount of mutagenic dNTPs [Zaccolo *et al.*].

Final PCR

For elimination of the mutagenic dNTPs, use an aliquot of 1 μ l of the first PCR reaction as template in a second PCR. Take 10x Mutagenesis

Buffer and add template, primers, dNTP Mix and Taq Polymerase in same concentrations as above. Fill up with PCR-grade Water to 50 μ l.

Use the same thermo cycling conditions but 20-30 cycles.

Selected References:

Zang *et al.* (2006) Efficient and high fidelity incorporation of dCTP opposite 7,8-dihydro-8-oxodeoxyguanosine by sulfolobus solfatarius DNA polymerase Dpo4. *Journal of Biological Chemistry* **281**:2358.

Zaccolo *et al.* (1999) The effect of high-frequency random mutagenesis on in vitro protein evolution: a study on TEM-1 beta-lactamase. *J. Mol. Biol.* **285 (2)**:775.

Crameri *et al.* (1998) DNA shuffling of a family of genes from diverse species accelerates directed evolution. *Nature* **391**:288.

Zaccolo *et al.* (1996) An approach to random mutagenesis of DNA using mixtures of triphosphate derivatives of nucleoside analogues. *J. Mol. Biol.* **255**:589.

Stemmer (1994) Rapid evolution of a protein in vitro by DNA shuffling. *Nature* **370**:389.

Cadwell et al. (1992) Randomization of genes by PCR mutagenesis. PCR Meth. Appl. 2:28.



Fig. 1: Structures of dPTP and 8-Oxo-dGTP





Fig. 2: Rate of mutagenesis as a function of the number of PCR-cycles [Zaccolo et al.]

Biorbyt Ltd. 5 Orwell Furlong, Cowley Road,Cambridge, Cambridgeshire CB4 0WY, United Kingdom Email: info@biorbyt.com | Phone: +44 (0)1223 859 353 | Fax: +44(0)1223 280 240 Biorbyt LLC Suite 103, 369 Pine Street, San Francisco California 94104, United States Email: info@biorbyt.com | Phone: +1 (415)-906-5211 | Fax: +1 (415) 651 8558